


**AN 434**

## Characterizing Annealed ULE B Implants using EAGLABS<sup>SM</sup> PCOR-SIMS<sup>SM</sup>

September 17, 2008 (Version 4.0)

### Introduction

Development and improvement of Ultra Low Energy (ULE) boron ion implantation is an area of intense interest as device dimensions continually shrink. Characterization of these implants requires accurate profile shape and oxide layer thickness within the upper several nanometers of the wafer surface. **EAGLABS<sup>SM</sup> PCOR-SIMS<sup>SM</sup>** represents the latest improvements in ULE B characterization that incorporates point-by-point data corrections for all regions of the profile. This method avoids near-surface profile distortions introduced by the older oxygen flooding and normal incidence techniques and yields the most accurate junction depth measurements due to precise measurement of surface oxide thickness.

### Discussion

The difference between old and new protocols is dramatically shown in the profiles above of a 250eV as-implanted and annealed Si wafer. Near surface artifacts in the old O<sub>2</sub>-flood and normal incidence protocols severely distort the B profile shape of both the as-implanted and annealed profiles. In contrast, the **EAGLABS<sup>SM</sup> PCOR-SIMS<sup>SM</sup>** protocol shows that the as-implanted sample has a Gaussian-shaped peak only 1.3nm below the surface. The annealed sample shows redistribution of B to the interface between the surface oxide layer and the Si substrate, in agreement with existing thermodynamic diffusion models at the Si/SiO<sub>2</sub> interface<sup>1</sup>. In addition, the **EAGLABS<sup>SM</sup> PCOR-SIMS<sup>SM</sup>** profile also gives a quantitative measure of the surface oxide thickness, a feature that is lacking entirely from the O<sub>2</sub>-flood and normal incidence profiles. Note that there has been little change in the oxide thickness due to annealing despite significant diffusion of B.

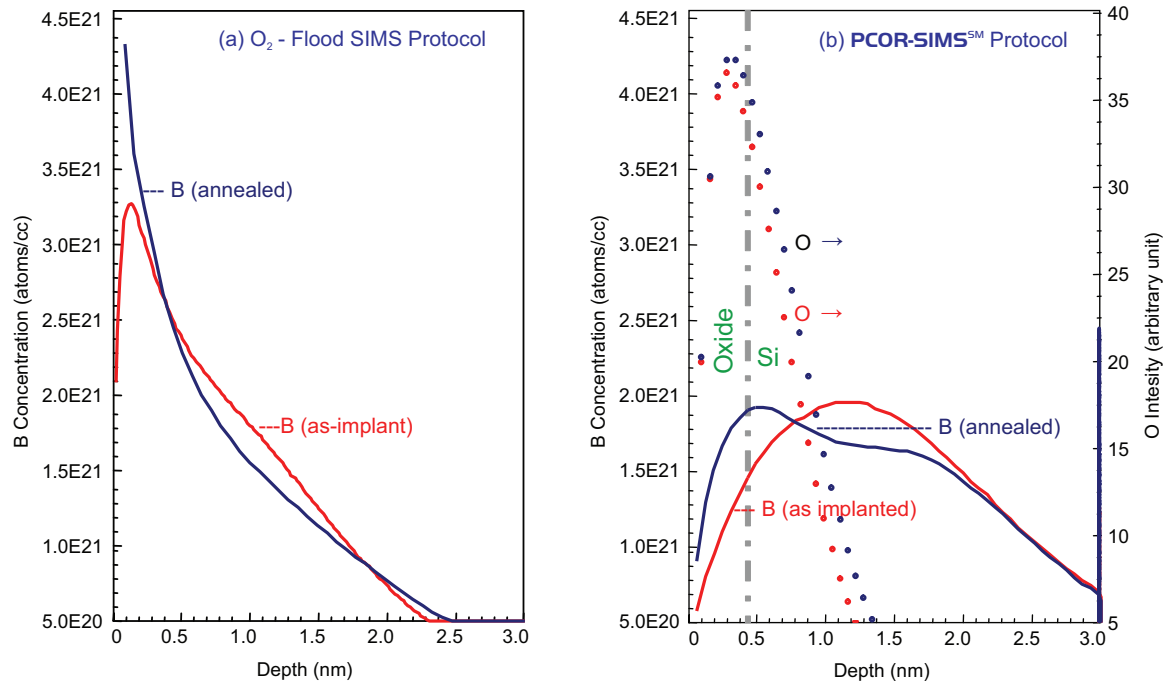


Figure 1. Comparison of two SIMS analysis protocols for 250eV boron implant characterization before and after anneal. Please note that concentration axes are linear unlike the usual log scales. **EAGLABS<sup>SM</sup> PCOR-SIMS<sup>SM</sup>** analysis of annealed B implantation detects accumulation at oxide/Si interface as expected (b). Whereas O<sub>2</sub>-flood SIMS protocol results in un-realistic profile shapes (a).

1 Phys. Rev. B 68, 195311 (2003)

\* The new PCOR-SIMS<sup>SM</sup> for ULE B protocol is the result of extensive development efforts by EAG. The "PCOR-SIMS<sup>SM</sup>" name describes, in part, EAG's proprietary methodology that includes point-to-point correction resulting in the most accurate SIMS profiling yet for ultra shallow implants.

---

Visit [www.eaglabs.com](http://www.eaglabs.com) for more information about all of EAG's services and solutions.

EAG Limited Administrative Offices, 810 Kifer Road, Sunnyvale, CA 94086 phone: 408 530 3500

Copyright © 2008 EAG Limited. All rights reserved. EAG, the EAG logo, are registered trademarks of EAG Limited.

Evans Analytical Group, Charles Evans & Associates, Thin Film Analysis, Inc., Applied Microanalysis Labs, Inc., AMIA Labs, Advanced Materials Engineering Research, Cascade Scientific Ltd., Cascade Scientific GmbH, Nano Science Corporation, Shiva Technologies, Inc., Shiva Technologies Europe SAS, Accurel Systems International Corporation, Micro Electronic Failure Analysis Services, Inc., DSL Labs Inc., White Mountain Labs LLC, are service marks of EAG Limited. All other company, product and service names may be trademarks of their respective companies. While every effort is made to ensure the information given is accurate, EAG Limited does not accept liability for any errors or mistakes which may arise. All information is subject to change without notice.